
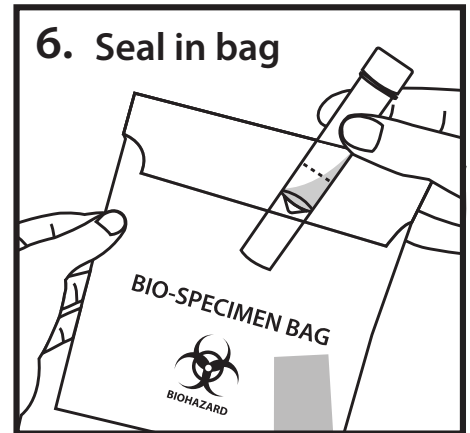
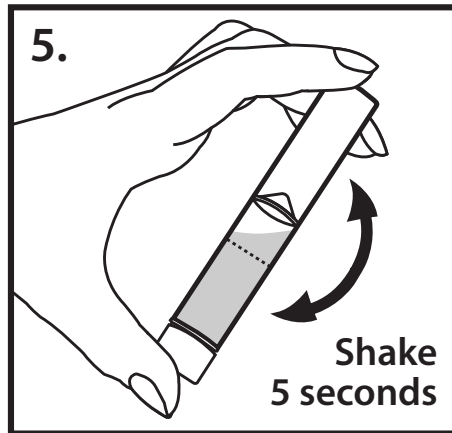
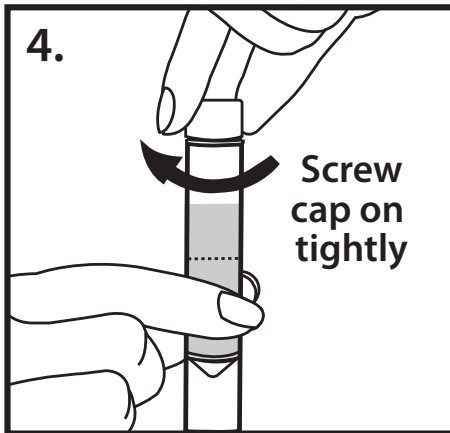
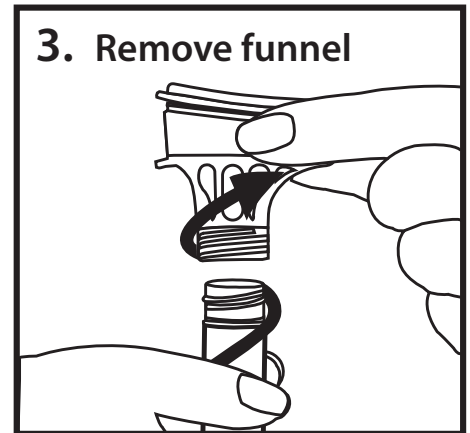


Oragene®-DISCOVER self-collection kit instructions

For complete instructions and additional languages, see inside kit.

Collection precautions: Do NOT eat, drink, smoke or chew gum for 30 minutes before giving your saliva sample.
Do NOT remove the plastic film from the funnel lid.

Procedure: Most people take between 2 and 5 minutes to deliver a saliva sample following steps 1 to 6.

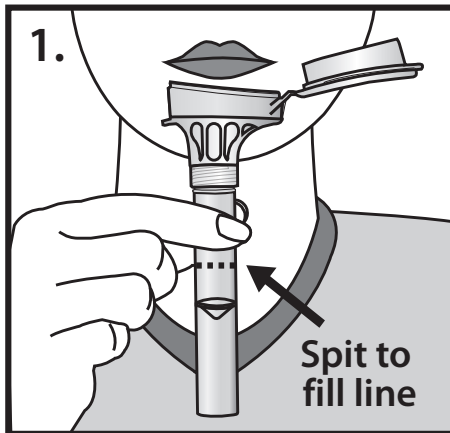

This product is designed for the collection of human DNA from saliva samples.

Warnings and precautions: Wash with water if the stabilizing liquid comes in contact with eyes or skin. Do NOT ingest.
See MSDS at www.dnagenotek.com.

Small cap, choking hazard.

Storage: Store at room temperature 15–30°C.

Oragene®-DISCOVER is for research use only, not for use in diagnostic procedures.
Some DNA Genotek products may not be available in all geographic regions.
*Oragene is a registered trademark of DNA Genotek Inc.