
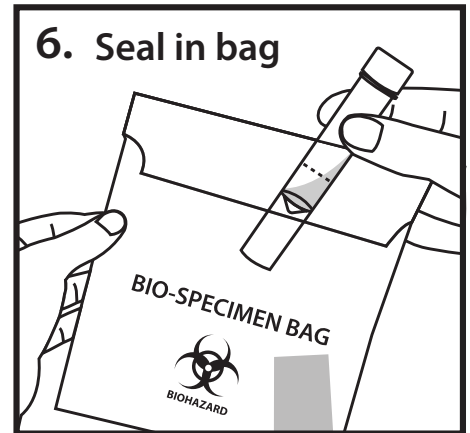
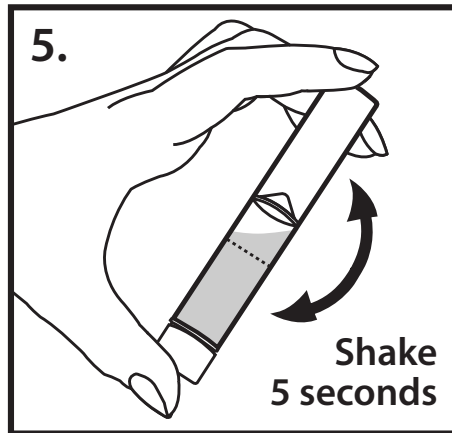
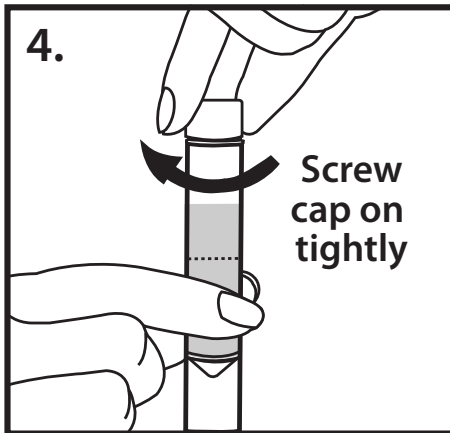
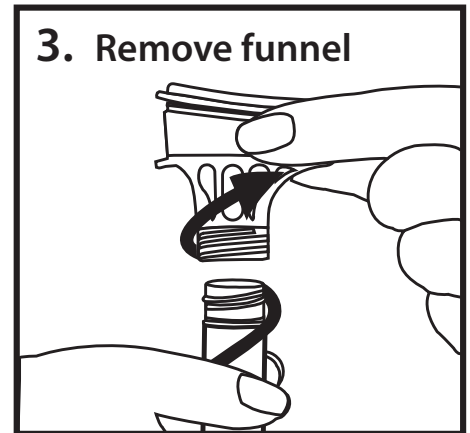


## Oragene®-DISCOVER self-collection kit instructions

**For complete instructions and additional languages, see inside kit.**

**Collection precautions:** Do NOT eat, drink, smoke or chew gum for 30 minutes before giving your saliva sample.  
Do NOT remove the plastic film from the funnel lid.

**Procedure:** Most people take between 2 and 5 minutes to deliver a saliva sample following steps 1 to 6.

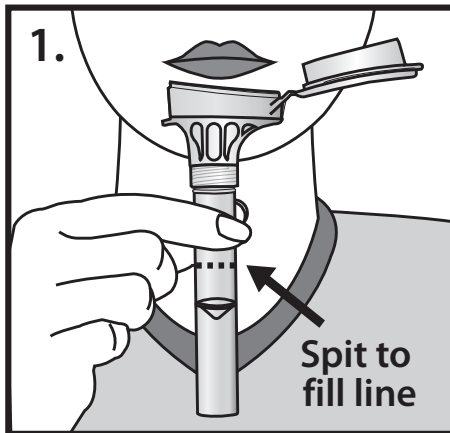

This product is designed for the collection of human DNA from saliva samples.

**Warnings and precautions:** Wash with water if the stabilizing liquid comes in contact with eyes or skin. Do NOT ingest.  
See MSDS at [www.dnagenotek.com](http://www.dnagenotek.com).

Small cap, choking hazard.

**Storage:** Store at room temperature 15–30°C.

Oragene®-DISCOVER is for research use only, not for use in diagnostic procedures.  
Some DNA Genotek products may not be available in all geographic regions.  
\*Oragene is a registered trademark of DNA Genotek Inc.