









Oragene®-DISCOVER self-collection kit instructions

For complete instructions and additional languages, see inside kit.

Collection precautions: Do NOT eat, drink, smoke or chew gum for 30 minutes before giving your saliva sample.

Do NOT remove the plastic film from the funnel lid.

Procedure: Most people take between 2 and 5 minutes to deliver a saliva sample following steps 1 to 6.

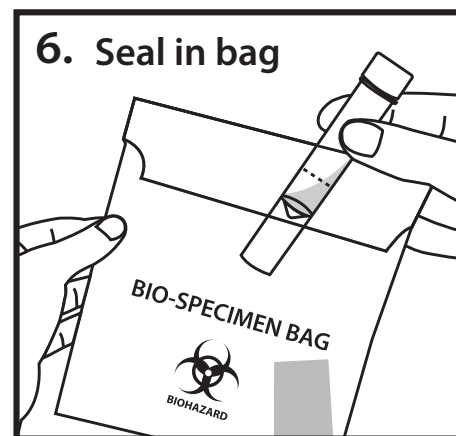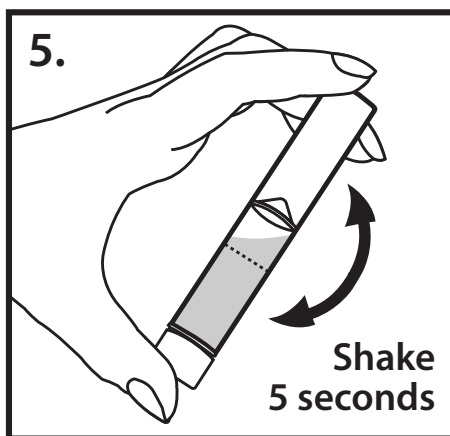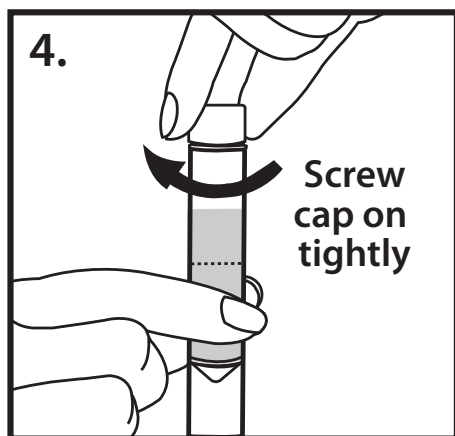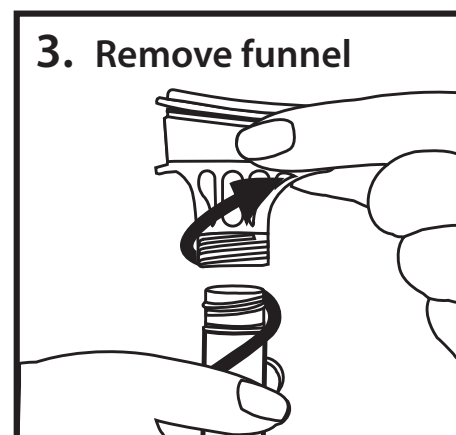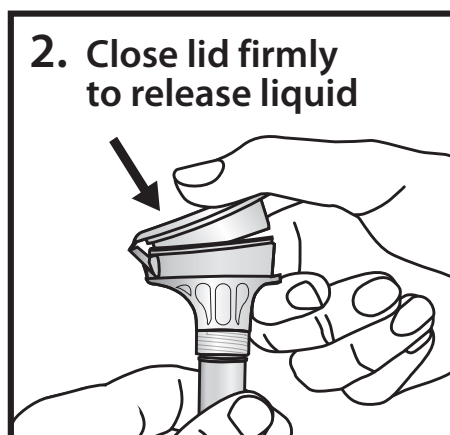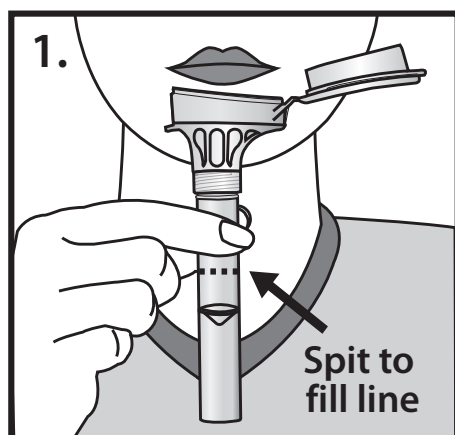

This product is designed for the collection of human DNA from saliva samples.

Warnings and precautions: Wash with water if the stabilizing liquid comes in contact with eyes or skin. Do NOT ingest. See MSDS at www.dnagenotek.com.

Small cap, choking hazard.

Storage: Store at room temperature 15–30°C.

Oragene®-DISCOVER is for research use only, not for use in diagnostic procedures. Some DNA Genotek products may not be available in all geographic regions. *Oragene is a registered trademark of DNA Genotek Inc.

ATSDR estimates the average public reporting burden for this collection of information as 10 minutes per response, including the time for reviewing instructions, searching existing data/information sources, gathering, and maintaining the data/information needed, and completing and reviewing the collection of information. An agency may not conduct or sponsor, and a person is not required to respond to a collection of information unless it displays a currently valid OMB Control Number. Send comments regarding this burden estimate or any other aspect of this collection of information, including suggestions for reducing this burden to CDC/ATSDR Information Collection Review Office, 1600 Clifton Road NE, MS D-74, Atlanta, Georgia 30333; ATTN: PRA (0923-0041).